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p94 LogTag Analyzer User Guide

## Manual USB driver installation on Windows XP

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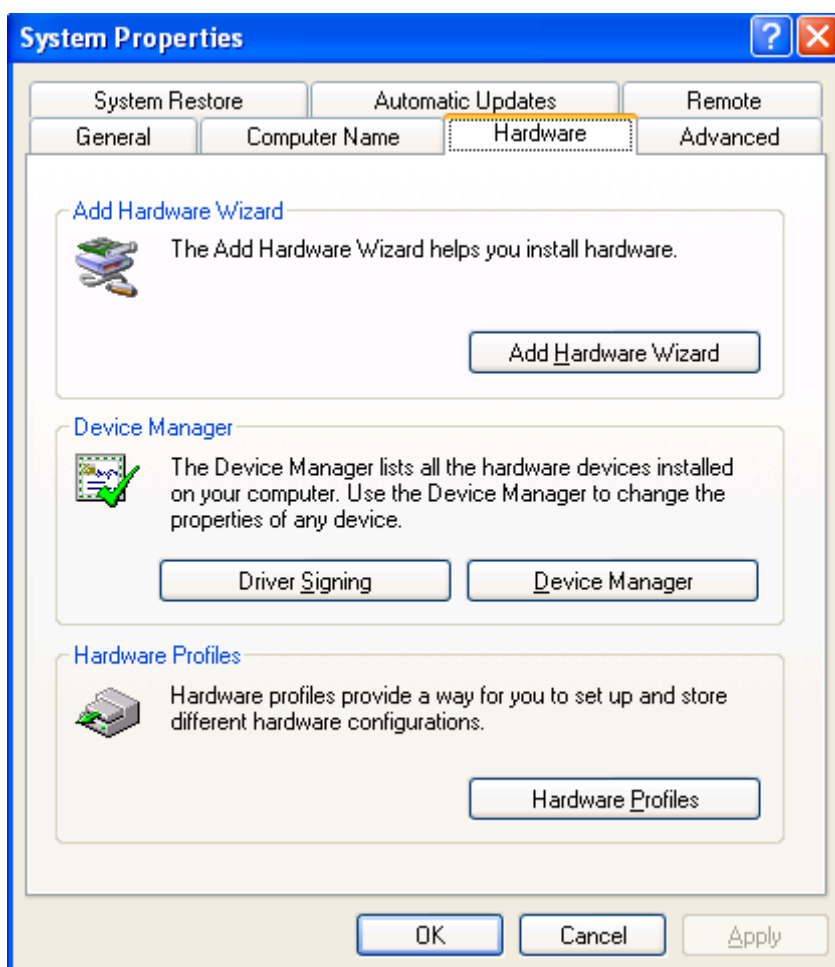
To manually reinstall the USB driver files, for Windows XP computers, open the "Device Manager". The "Device Manager" can be started by opening the "System" icon, located in

the Windows "Control Panel", selecting the "Hardware" tab in the "System Properties" window and then click the "Device Manager" button.


## Chapter 11 Appendix p95

1 Using the mouse (pointing device), click the right button to display the popup context menu, which will be similar to the following picture.

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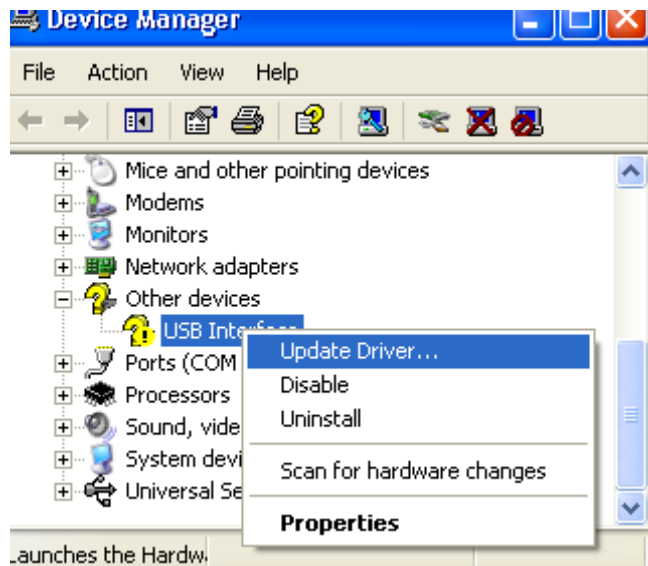2 Select the "Update Driver..." option, using the left button on the mouse, which will cause the device driver wizard to start, similar to the following picture. Ensure the "Install from a list or specific location (Advanced)" option and then click the "Next >" button to continue.

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## p96 LogTag Analyzer User Guide

3 When prompted for the location of the driver files, enter the location of the installation of the software followed by the path to the USB driver files. The default location for these driver files is "C:\Program Files\LogTag Recorders\LogTag Analyser\Drivers\USB Interface 2.0". You may use the "Browse..." button in the window that appears, similar to the following picture to help you locate the appropriate folder.

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
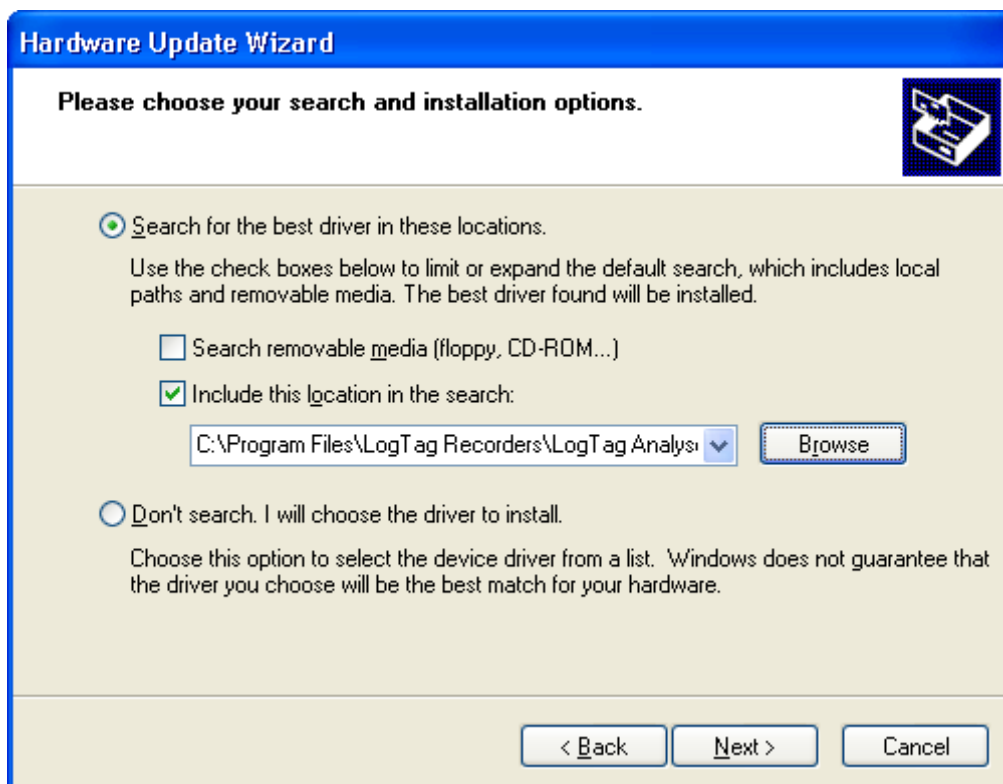
## Chapter 11 Appendix p97

4 Once the driver files have been installed successfully you will see a window similar to the following picture. You may now connect and start using your USB Interface Cradle(s).

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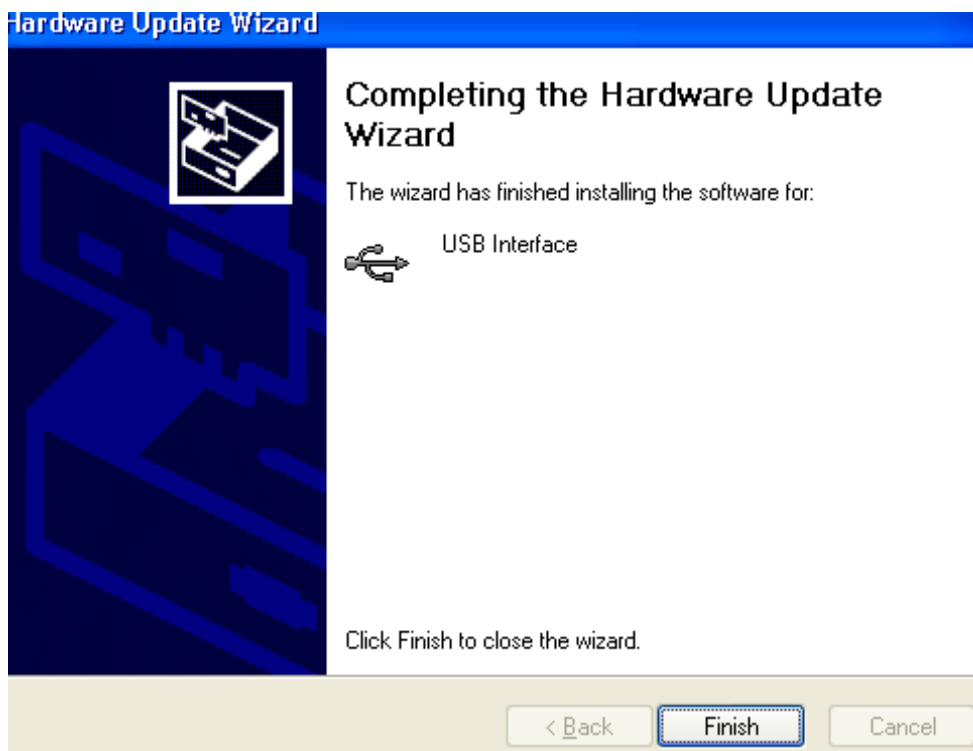Note, if you see a window similar to the following picture appear during the installation of the USB driver files, you must click the "Continue Anyway" button for the USB driver files to successfully install.

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The screenshot shows the LogTag configuration window with the following settings:

- Action List:**
  - 1. Welcome and general instructions
  - 2. Locate LogTag(s), download and save relevant data
  - 3. Configure LogTag(s) for next use
  - 4. Prepare LogTag(s) for next use
- User ID:** Description of monitored environment
- Start Method:** Push button start
- Logging Options:**
  - Enable pre-start logging
  - Record readings continuously, overwrite oldest when memory full
  - Record readings so that:
    - Readings recorded will span at least 25 days
    - Number of readings to record 7,200 (maximum is 8003)
- Recording Interval:** Record a reading every 5 Minutes
- Delay:** Begin recording after a delay of 0 Minutes
- Enable the OK (Green) indicator

**Alert Settings (Humidity/Temperature):**

- Enable the Alert (Red) indicator, when:
  - Readings below 32.6 or above 58.2 %RH.
  - After 24 consecutive alert readings (2 Hours)
  - After 0 alert readings have occurred (None)
  - Leave alert indicator enabled even if readings return within alert range
- Clear and reset alert when START MARK button pressed
- Configure requires a password

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